

Metallographic preparation of fasteners




Struers

Application Notes

Pins and rods as means of mechanically connecting parts are one of the oldest construction elements. Rivets and screws were first made from precious metals and later from non-ferrous metals and steels through forging and mechanical cutting. Not until the development of metals that were ductile enough to withstand the pressure of pressing and rolling tools, was cold forming adopted as a process used for the mass production of fasteners.

Both, rivets and screws have their specific fields of applications, such as aircraft and automotive industry, that are determined by specifications, safety and economical considerations.

In many engineering applications, fasteners play a crucial role with regard to structural safety, and therefore have to be manufactured in compliance with specific standards. As a consequence, process and quality control procedures in general and metallographic examination in particular, are important in the production process. The tests range from checking dimensions, mechanical properties, and physical variables, to structure examination and quantitative analysis. Metallographic examination is part of the quality control regimen for producing fasteners, may it be spot checks of raw material, control and verification of forming and heat treatment processes or failure analysis.

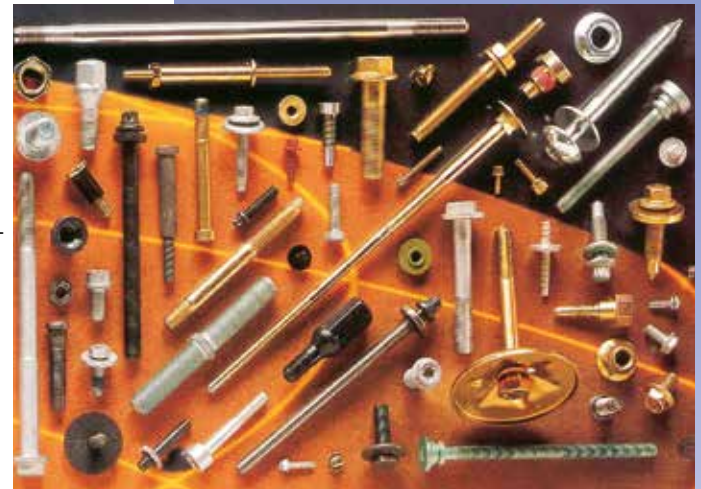

Riveted joint, steel rivet in aluminium

The main categories of mechanical joining are threaded fasteners such as bolts, nuts and screws, and non-threaded fasteners, like rivets and pins. Depending on the requirements regarding strength, torsion, environmental conditions etc. threaded and non-threaded fasteners are used for different applications.

Difficulties during metallographic preparation:


Cutting: the size and shape of fasteners does not always allow for secure clamping and proper cross sectional cutting.

Mounting: shrinkage gaps on the thread and head of a fastener can make it difficult to achieve good edge retention and clean the sample properly in order to observe it microscopically.


Solution:

Cutting: Special holder for fasteners, precision cutting for small screws, or mounting of complete fastener with subsequent coarse grinding to the center.

Mounting: Proper degreasing and cleaning, use of mounting media material with the least shrinkage for hot compression or cold mounting.

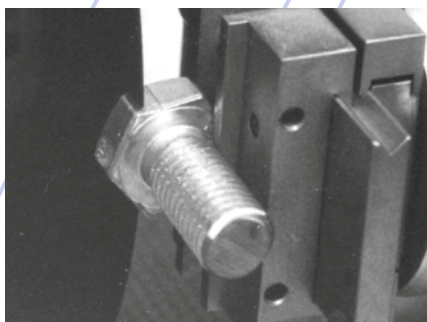

Cutting excess of screw head off prior to mounting

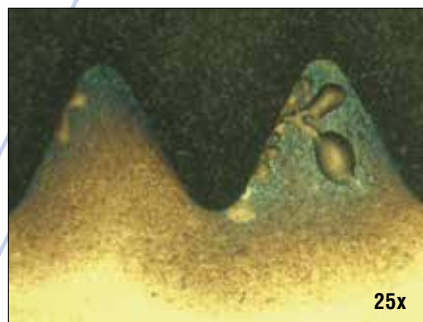

Thread, staining due to shrinkage gap

Production and application of fasteners

In the historical development of fastener production mechanical machining was important, especially for cutting threads. Today, chipless shaping (rolling) is a universal and economical production method, and machining is only used after the initial shaping, for finishing very special, high quality construction screws which require specific shapes, tolerances and surfaces.


Forming is carried out through either cold or hot pressing. Cold forming technology and metallurgy have advanced to a state where cold pressing has almost replaced hot forming in the production of fasteners.

For very large or complex shaped bolts, hot forming is still an option. Large section sizes, necessitate high cold forming forces resulting in an increased severity of the effects of cold work. For cold forming a continuous wire is fed into an extrusion press, where through compression and reduction of the wire diameter in a die the fastener is shaped (Fig. 1). The threads are formed through a cold rolling operation (Fig. 2). As the threads are the most crucial part on a screw, correct rolling operation and subsequent heat treatment are very important. The cold working processes increase the hardness, and for quality steel fasteners, various heat treatment steps are carried out to achieve certain material properties for specific applications. Low carbon steels are carburized so that the centre is soft and ductile and the surface hard. Carbon steels are hardened and tempered, and depending on the application, temperature variations for austenitization and tempering are used to achieve different mechanical properties from the same raw material. Additional surface hardening, for instance on the heads, can be achieved through induction hardening. During the various stages of shaping and heat treating other steps such as pickling, cleaning and oiling take place.

Coating for corrosion protection is a last and usually a separate step from the actual manufacturing. For improved corrosion protection, fasteners are coated with manganese, zinc or iron phosphate and then oiled.

Fig. 1. Cold Heading.


A basic operation, it forms round wire stock into the essential various shapes while still "cold". A continuous, automatic operation it nonetheless must retain continuity of grain to assure the full integrity of the fastener.

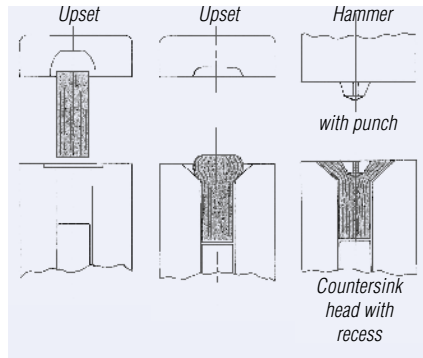

Galvanizing with zinc and cadmium or chromium plating offer a higher corrosion protection, and so do vapor deposition coatings of cadmium and aluminium. (In some countries cadmium coatings are no longer applied for environmental reasons.)

In addition to low carbon and carbon steel fasteners for general applications, fasteners are made of high strength steel for construction and mechanical engineering applications, stainless steel for corrosion resistance, nickel and cobalt based alloys for high temperature applications, and titanium in the aircraft, medical and food industry. Non ferrous fasteners of brass and copper are used in connection with decorative metal rails and trimmings. Steel rivets are used in the automotive industry and titanium rivets are mainly used in aircraft outer body construction .

A main prerequisite for manufacturing good fasteners is quality raw material. There can be defects in the bars such as seams, voids in the centre of the bar, or incorrect heat treatment that produces the wrong grain size or structure. These can all lead to defects during the production of the fasteners. Metallographic quality control of raw material is therefore as important as the general production control.

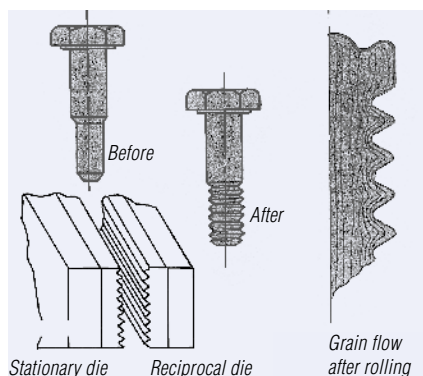

Difficulties in the preparation of fasteners

The challenge of metallographic preparation of fasteners is their geometry. Usually the cut has to go through the centre of the screw and as the head is a protruding part, it can be difficult to clamp a rivet or screw securely to make a cross sectional cut. Large bolts can usually be cut in half without difficulties, however the thinner the screw is, the more difficult it is to cut. The problem with geometry also influences the mounting, as the curves of thread and head are spots where shrinkage of mounting resin can occur preferentially. This is particular crucial on coated material, because coatings can not be examined properly without good edge retention.



Recommendations for the preparation of fasteners

To overcome the problems of cutting and mounting the following can be recommended:

For medium sized or thin screws a special holder with threads can be made that holds the part securely during cross sectioning (Fig. 3). For smaller screws or rivets the head can be cut off on one side, so that the screw can be laid flat into the mounting press. After mounting, the rest of the screw can be ground to the centre. An alternative is a special lower ram for the mounting press that has a recess on both sides into which the screwheads fit (Fig. 4). For very small screws cutting is not efficient. They should be mounted with a piece of metal or plastic as a shim to level the screw and then be ground to the middle after mounting (Fig. 5).



Fig. 3: Special holder for cutting screws

Fig. 2: Thread rolling.

A blank is rolled between two flat dies with precise parallel grooves of the thread type required. A cold - or warm - forming operation, rolled threads are strong and smooth - and no material is wasted.

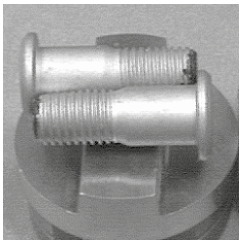


Fig. 4

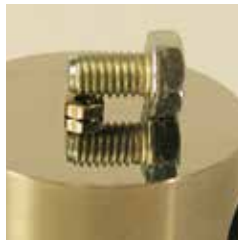


Fig. 5

Fig. 4:
Special lower ram for hot
compression mounting

Fig. 5:
Shimming of screw for
mounting

The most efficient way of grinding fasteners is with a grinding stone on an automatic grinding machine, especially if large volumes have to be handled.

Dependent of material properties, for hot compression mounting or resins containing harder filler material (DuroFast) phenolic resin with carbon filler (PolyFast) or melamine wite mineral and glass filler (LevoFast) are recommended. For cold mounting epoxy resins (Epofix) epoxies are recommended as they have the least shrinkage. Before mounting, the parts should be thoroughly cleaned, with a degreasing agent such as acetone or ethanol. Particular attention should be paid to the threads and the radius below the head to make sure that the resin will adhere properly to the sample material.

Grinding and polishing

Once the fasteners are properly cut and mounted, the grinding and polishing has to be carried out according to the specific material. The wide range of fastener materials can not all be covered in this Application Note. We have selected two types of metal fasteners of which the preparation methods are given in tables 1 and 2. (For more detailed information on preparations, please refer to the Struers e-Metalog or Struers Application Notes). The examples for preparation data given in this Application Note are for brass shrews (Table 1) and screws of low alloyed steel (Table 2). The data are for 6 mounted samples, 30 mm dia., clamped into a holder. The coarse grinding to the middle of the screws was carried out on the Struers automatic grinding machine Abraplan, the subsequent automatic fine grinding and polishing on Struers Tegramin.



Etching and structure interpretation

For etching of metallographic sections of fasteners the common, recommended etching solutions for the respective materials, such as carbon steel, titanium, brass, stainless steel etc. are used.



Fig. 6:
Fatigue fracture

As mentioned above, during the production of fasteners mechanical defects such as burrs, cracks (Fig. 7a), folds and overlaps can occur, some of which can be detected through visual inspection or with non-destructive methods such as dye penetrant and magnetic particle inspection. The depth of these defects can be detected by metallographic means (Fig. 7b). Microstructural defects are mostly due to heat treatment such as overheating (Fig. 8), grain growth, intergranular oxidation and decarburisation (Fig. 9).

In many applications, the fracture of threaded fasteners can have serious consequences, consequently metallographic structural analysis is essential to find out what caused the fracture. They can occur through certain influences over time, such as mechanical, like bending and straining, or thermal and environmental exposure to lubricants, steam, chemical reagents etc. This can lead to fractures due to fatigue (Fig.6), torsion, embrittlement, corrosion and other multiple causes.

Summary

The term "fasteners" includes a wide range of threaded and non threaded bolts, screws, rivets and pins which all require specific raw material



Fig. 7a: Split in head of carbon steel bolt caused by a seam.



Fig. 7b: Traverse section of the seam in the fig. 7a. Echant: 3% Nital

Preparation Method

Brass screw

Grinding


Step	PG	FG	
Surface	Foil/Paper	MD-Largo	
Abrasive	Type	SiC	Diamond
	Size	#320	9 µm
Suspension/Lubricant	Water	DiaPro Allegro/Largo 9	
rpm	300	150	
Force [N]/specimen	30*	30*	
Time (min)	As needed	3	

Polishing

Step	DP	OP	
Surface	MD-Mol	MD-Chem	
Abrasive	Type	Diamond	Collodial Silica
	Size	3 µm	0.04 µm
Suspension/Lubricant	DiaPro Mol R 3	OP-S NonDry**	
rpm	15	150	
Force [N]/specimen	30*	15*	
Time (min)	3	3	

Table 1

* Value for 6 mounted samples, 30 mm diam. Clamped in a holder.
**96 ml OP-S NonDry
2 ml ammonia (25%)
2 ml hydrogen peroxide (33%)

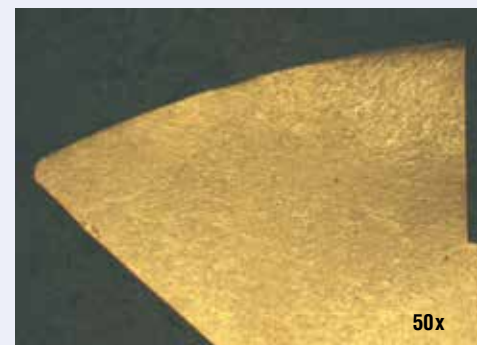

Head of brass screw with flow lines from cold forming. Echant: Iron (III) chloride

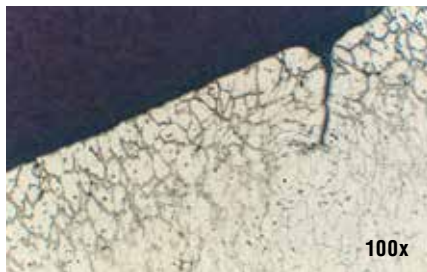


Fig. 8:
Melted grain boundaries
due to overheating in heading.
Nickel base alloy
Etchant:
modified Kalling's

Struers ApS
Pederstrupvej 84
DK-2750 Ballerup, Denmark
Phone +45 44 600 800
Fax +45 44 600 801
struers@struers.dk
www.struers.com

Preparation Method


Bolt, alloyed steel
Grinding

Step		PG	FG
	Surface	Stone	MD-Allegro
	Abrasive	Type	Aluminium oxide
		Size	#150
	Suspension/ Lubricant	Water	DiaPro Allegro/Largo 9
	rpm	1500	150
	Force [N]/ specimen	35*	35*
	Time (min)	As needed	4

Polishing

Step		DP	OP
	Surface	MD-Dac	MD-Nap
	Abrasive	Type	Diamond
		Size	3 µm
	Suspension/ Lubricant	DiaPro Dac 3	DiaPro Nap B 1
	rpm	150	150
	Force [N]/ specimen	35*	25*
	Time (min)	4	1-5

Table 2 * Value for 6 mounted samples,
30 mm in diam. clamped in
a holder.

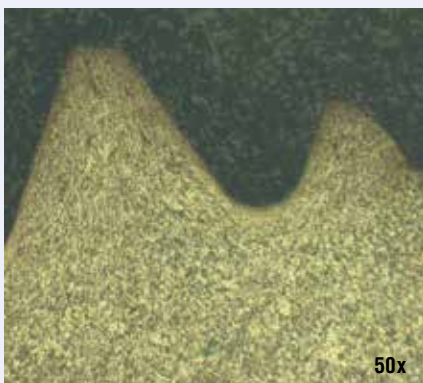

Thread of bolt, alloyed steel tempered
Etchant: 3% Nital

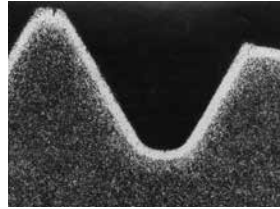


Fig. 9:
Section of thread
showing
decarburization
Etchant:
5% Nital

and production processes. Good metallographic quality control for raw material and production is essential to ensure quality fasteners. The metallographic checks are mainly for mechanical and thermal damages due to production and subsequent heat treatment. Interpretation of structure and defects requires some experience as the range of materials and shapes of fasteners is very wide.

Specific problems during the metallographic preparation are cutting and mounting, which can be overcome by using various helpful tools. The grinding and polishing procedure depends on the fastener material to be prepared and can be carried out on automatic equipment with a three or four step method.



Application Notes

Metallographic preparation of fasteners
Elisabeth Weidmann, Struers A/S, Denmark
John McElwain, Struers Inc., Westlake OH, USA

Acknowledgement

We wish to thank Frauke Hogue, Hogue Metallography, Pacific Palisades, Ca, USA, for contributing information on microstructures of fasteners and the permission to reproduce figs no. 7a+b and 8.

Figs. 1 and 2: Copyright 1975 Alcoa Global Fasteners, Inc. and its affiliated companies. All rights reserved.

Illustrations are re-printed by permission from copyright owner. Reproduction of the illustrations in whole or in part without written permission of the copyright owner is strictly prohibited.

Bibliography

Handbuch der hochfesten Schrauben
K.-H. Kübler, W. J. Mage, Hrsg.: Kamax-Werke,
Verlag W. Giradet, Essen, 1986

Schrauben Vademecum, K. H Illgner, D. Blume,
Bauer und Schauer, Neuss, 1978

Characterization of defects in precision fasteners by
metallographic methods.

Frauke Hogue, Structure 30, 1996

AUSTRALIAN & NEW ZEALAND
Struers Australia
27 Mayneview Street
Milton QLD 4064
Australia
Phone +61 7 3512 9600
Fax +61 7 3369 8200
info.au@struers.dk

BELGIUM (Wallonie)
Struers S.A.S.
370, rue du Marché Rollay
F-94507 Champigny
sur Marne Cedex
Téléphone +33 1 5509 1430
Télécopie +33 1 5509 1449
struers@struers.fr

BELGIUM (Flanders)
Struers GmbH Nederland
Zomerdijk 34 A
3143 CT Maassluis
Telefoon +31 (10) 599 7209
Fax +31 (10) 5997201
netherlands@struers.de

CANADA
Struers Ltd.
7275 West Credit Avenue
Mississauga, Ontario L5N 5M9
Phone +1 905-814-8855
Fax +1 905-814-1440
info@struers.com

CHINA
Struers Ltd.
No. 1696 Zhang Heng Road
Zhang Jiang Hi-Tech Park
Shanghai 201203, P.R. China
Phone +86 (21) 6035 3900
Fax +86 (21) 6035 3999
struers@struers.cn

CZECH REPUBLIC & SLOVAKIA
Struers GmbH Organiza ni slozka
v dektechnicky park
P ílepská 1920,
CZ-252 63 Roztoky u Prahy
Phone +420 233 312 625
Fax +420 233 312 640
czechrepublic@struers.de
slovakia@struers.de

GERMANY
Struers GmbH
Carl-Friedrich-Benz-Straße 5
D- 47877 Willich
Telefon +49 (0) 2154 486-0
Fax +49 (0) 2154 486-222
verkauf@struers.de

FRANCE
Struers S.A.S.
370, rue du Marché Rollay
F-94507 Champigny
sur Marne Cedex
Téléphone +33 1 5509 1430
Télécopie +33 1 5509 1449
struers@struers.fr

HUNGARY
Struers GmbH
Magyarország Fióktelep
Tatai út 53
2821 Gyermely
Phone +36 (34) 880546
Fax +36 (34) 880547
hungary@struers.de

IRELAND
Struers Ltd.
Unit 11 Evolution@ AMP
Whittle Way, Catcliffe
Rotherham S60 5BL
Tel. +44 0845 604 6664
Fax +44 0845 604 6651
info@struers.co.uk

ITALY
Struers Italia
Via Monte Grappa 80/4
20020 Arese (MI)
Tel. +39-02/38236281
Fax +39-02/38236274
struers.it@struers.it

JAPAN
Marumoto Struers K.K.
Takanawa Muse Bldg. 1F
3-14-13 Higashi-Gotanda,
Shinagawa
Tokyo
141-0022 Japan
Phone +81 3 5488 6207
Fax +81 3 5488 6237
struers@struers.co.jp
13.04.2018 R04 / 62140203

NETHERLANDS
Struers GmbH Nederland
Zomerdijk 34 A
3143 CT Maassluis
Telefoon +31 (10) 599 7209
Fax +31 (10) 5997201
netherlands@struers.de

NORWAY
Struers ApS, Norge
Sjøskogenveien 44C
1407 Vinterbro
Telefon +47 970 94 285
info@struers.no

AUSTRIA
Struers GmbH
Zweigniederlassung
Österreich
Betriebsgebiet Puch Nord 8
5412 Puch
Telefon +43 6245 70567
Fax +43 6245 70567-78
austria@struers.de

POLAND
Struers Sp. z o.o.
Oddział w Polsce
ul. Jasnogórska 44
31-358 Kraków
Phone +48 12 661 20 60
Fax +48 12 626 01 46
poland@struers.de

ROMANIA
Struers GmbH, Sucursala Bucuresti
Str. Preciziei nr. 6R
062203 sector 6, Bucuresti
Phone +40 (31) 101 9548
Fax +40 (31) 101 9549
romania@struers.de

SWITZERLAND
Struers GmbH
Zweigniederlassung Schweiz
Weissenbrunnenstraße 41
CH-8903 Birmensdorf
Telefon +41 44 777 63 07
Fax +41 44 777 63 09
switzerland@struers.de

SINGAPORE
Struers Singapore
627A Aljunied Road,
#07-08 BizTech Centre
Singapore 389842
Phone +65 6299 2268
Fax +65 6299 2661
struers.sg@struers.dk

SPAIN
Struers España
Camino Cerro de los
Gamos 1
Building 1 - Pozuelo de
Alarcón
CP 28224 Madrid
Teléfono +34 917 901 204
Fax +34 917 901 112
struers.es@struers.es

FINLAND
Struers ApS, Suomi
Hietalahdenranta 13
00180 Helsinki
Puhelin +358 (0)207 919
430
Faksi +358 (0)207 919 431
finland@struers.fi

SWEDEN
Struers Sverige
Box 20038
161 02 Bromma
Telefon +46 (0)8 447 53 90
Telefax +46 (0)8 447 53 99
info@struers.se

UNITED KINGDOM
Struers Ltd.
Unit 11 Evolution @ AMP
Whittle Way, Catcliffe
Rotherham S60 5BL
Tel. +44 0845 604 6664
Fax +44 0845 604 6651
info@struers.co.uk

USA
Struers Inc.
24766 Detroit Road
Westlake, OH 44145-1598
Phone +1 440 871 0071
Fax +1 440 871 8188
info@struers.com